



User Guide


Iron Test Kit

Always remember

1. To wash the test vials properly before test.
2. While matching the colour of the test sample with colour, always place the vial against a white background, to get maximum accuracy.
3. Test the samples as soon as possible after collecting it from the source.

Test Procedure:

1. Fill the Test vial with 4 ml of pond/ tank water using syringe.
2. Add 4 drops of Iron Reagent 1, close the cap and shake for 5 seconds, followed by adding 4 drops of Iron Reagent 2 and shaking for 5 seconds. Finally add 4 drops of Iron Reagent 3.
3. Close the cap and mix rigorously for five seconds.
4. Leave the tube for two minutes and compare the colour with colour card.

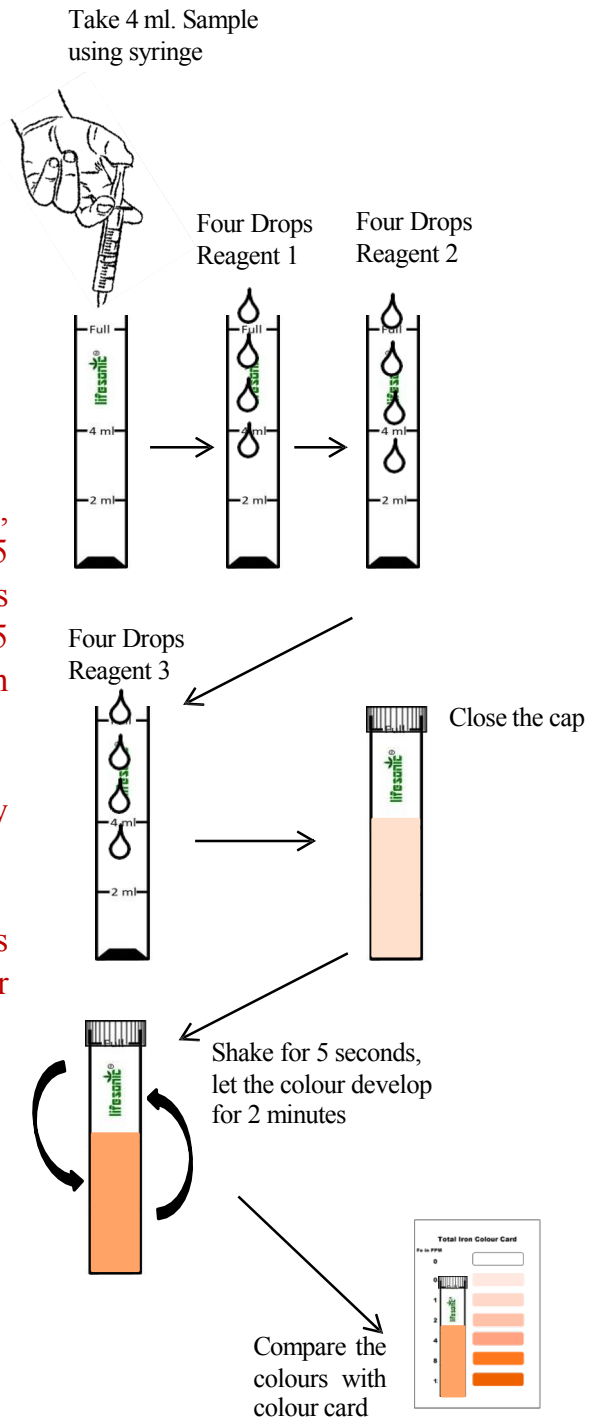

Store at 15-20°C in dry and dark place